

BioFX™ Alkaline Phosphatase/Horseradish Peroxidase Block

Product insert

Product number:

APHP-0100-01 (100 mL) Custom Sizes Available

Product use:

BioFX™ Alkaline Phosphatase/Horseradish Peroxidase Block is supplied as a ready-to-use blocking reagent for immunohistochemistry samples. The solution blocks both alkaline phosphatase and peroxidase endogenous enzymatic activity.

Product stability and storage:

Product stability	3 years from the date of manufacture.
Storage	Store at 2-8°C. Care should be taken to minimize the risk of contamination.

Recommendations for use:

- 1) Samples should be blocked for a minimum of 10 minutes at room temperature (25°C).
 - Additional blocking time may be required depending on tissue thickness and sample permeability.
- 2) Samples should be rinsed with copious amounts of distilled water to remove excess blocking solution.

For technical assistance, email ivdtechsupport@surmodics.com

Related Products:

BioFX® DAB Immunohistochemistry Substrate (DABI)

SURMODICS and BIOFX are trademarks of Surmodics, Inc. and/or its affiliates.

These products are for research and/or further manufacture and are not intended for use in humans or therapeutic purposes. Sales are without any seller's warranty or representation, expressed or implied, by usage or otherwise; no claims beyond replacement of unacceptable material or refund of purchase price shall be allowed. All claims must be made within 30 days following date of delivery.